

How to Customize settings in Nikon Scan

The Nikon Scan program has several settings that can be customized, but these must be manually saved.

How to change the default format for saved files:

- 1) While in the Nikon Scan scanning window, click on the blue button for “Preferences.” Choose “File Saving” from the left hand column and then select the desired option.
- 2) Click on “OK.” The setting should now be saved on your profile.

How to save options not listed under “Preferences” (i.e. filters, etc.)

Please note: Settings that are not listed under “Preferences” must be loaded each time you use the Nikon Scan program. Instructions for loading these are listed below

1. While in Nikon Scan, change any settings you need to customize.
2. Click on **Settings** and then **Export Settings** from the scanning window. Enter a filename and, if desired, specify a location. Click on **Save**.
3. You may exit Nikon Scan normally. If you receive a message stating “Unable to save settings file,” you may ignore it.

How to retrieve a saved settings file

1. While in the Nikon Scan scanning window, click on **Settings** and then **Import Settings**.
2. Browse to the location of your settings file and select it. Click on **Open**. Your settings should now be loaded.

If you have any problems saving settings in Nikon Scan, please see the lab assistant on duty for more help.